

How to upgrade the TrueChrome HDMI camera firmware

Firmware Upgrade steps:


1. Connect the SD card to a PC and create a folder “update” in the SD card root directory. Copy the firmware file “*.bin” to this “update” folder.



Note: The created folder name in the SD card ROOTDirectory MUSTbe “update”.

2. Insert the SD card into the camera SD card slot. Plug in the camera power, attach the USB mouse to the camera. Attach the camera to a monitor via HDMI cable and switch the monitor to HDMI input mode.

<Note: the monitor setting interface might be different because of different brand monitors.>

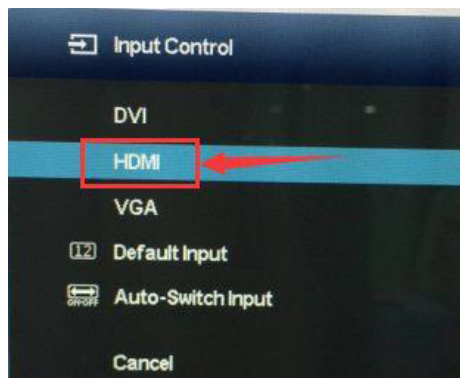

3. Switch on the camera power. A message “Detects new firmware Update” will be popped up in the preview screen, click on “Yes” to start the firmware upgrade.

Note: During the upgrade, please **DO NOT** power off the camera.

4. After finish the upgrade, a message “Update success restart now” shows up. Wait for the camera automatically restarting. Then if it goes to the preview images, it means the firmware is upgraded successfully.

