

MA502  
or  
MA504  
Eyepieces

MA555/05  
Gem Clamp

MA557  
Iris Diaphragm

GEMZ/PT  
Gemological Stand

EMZ-5TR  
Microscope Head

EMZ = zoom body

EMT = turret body



MA309/100

consists of

MA309/25 transformer  
and

MA309/PT#1 light fixture

GEMZ/PT = for zoom bodies only

GEMT/PT = for turret bodies only

## **Setting Up the GEMZ Stand**

Make sure all the plugs are plugged properly under the stage.

After plugging the unit:

The top light has a start-up button. When clicked once it acts as a starter (one bulb goes on). The second click gives the user a brighter intensity on the incident light (Black Box).

When the white on/off switch under the stage is turned off while the light is at high intensity, turning the switch back on will not turn the light on at all.

When the switch is turned off while the light is at low intensity, turning the switch back on will turn the light on at low intensity.

The silver knob on the bottom of the stand located on the left rear-off/on switch controls only the transmitted light portion of the stand. Make sure this is plugged properly.