



MEIJI TECHNO AMERICA

5895 Rue Ferrari, San Jose, CA 95138, USA


Tel : 408-226-3454 Fax : 408-226-0900

Toll Free : 1-800-832-0060

E-mail : info@meijitechno.com

http://www.meijitechno.com

Installation Diagram

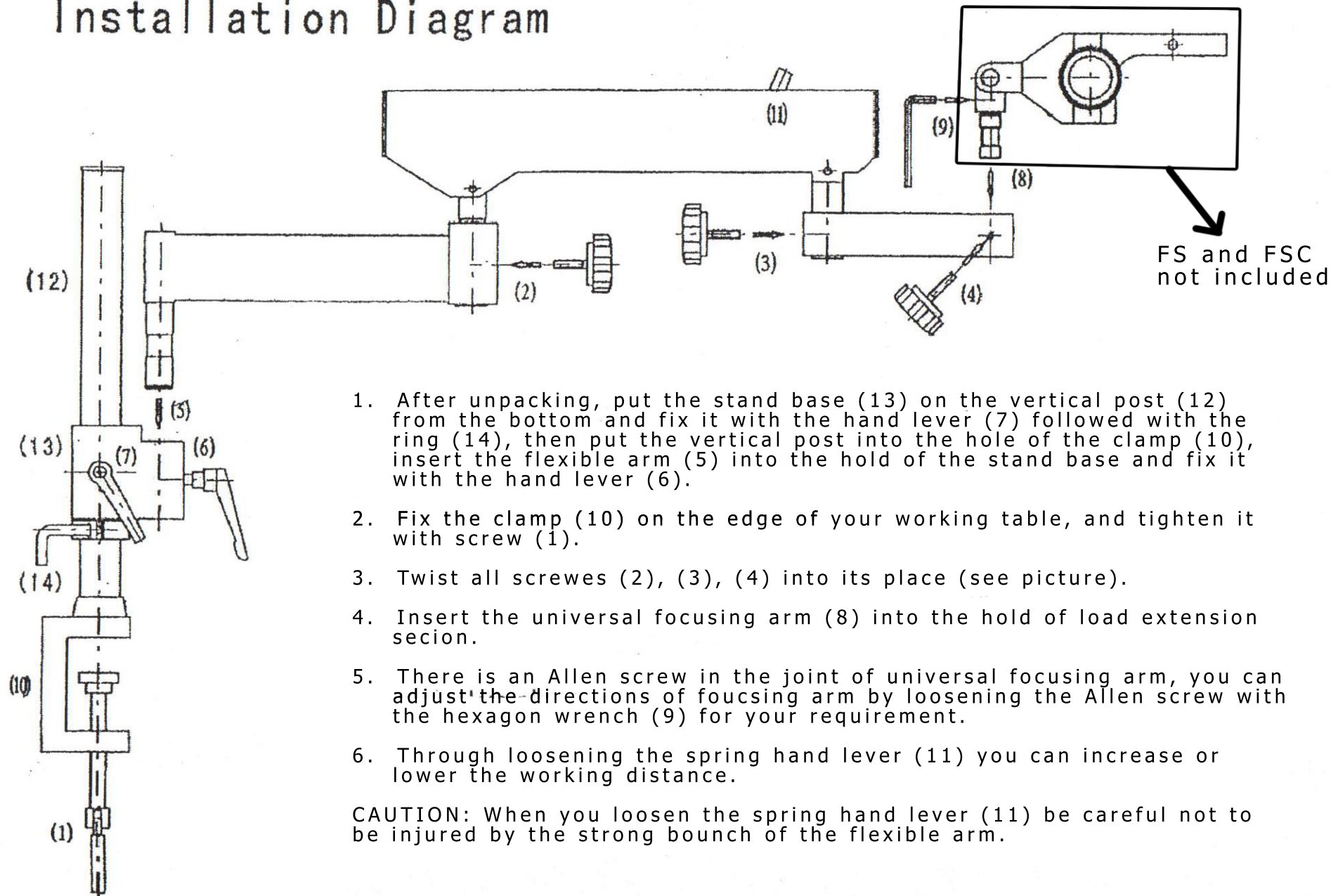

1. After unpacking, put the stand base (13) on the vertical post (12) from the bottom and fix it with the hand lever (7) followed with the ring (14), then put the vertical post into the hole of the clamp (10), insert the flexible arm (5) into the hold of the stand base and fix it with the hand lever (6).
2. Fix the clamp (10) on the edge of your working table, and tighten it with screw (1).
3. Twist all screws (2), (3), (4) into its place (see picture).
4. Insert the universal focusing arm (8) into the hold of load extension section.
5. There is an Allen screw in the joint of universal focusing arm, you can adjust the directions of focusing arm by loosening the Allen screw with the hexagon wrench (9) for your requirement.
6. Through loosening the spring hand lever (11) you can increase or lower the working distance.

CAUTION: When you loosen the spring hand lever (11) be careful not to be injured by the strong bounce of the flexible arm.