

Micro WiFi EDU Troubleshoot Instructions

1. Switch off all the related components and software, Apps
2. The first step is to switch off the Master Router, unplug it. Wait for 20 seconds, and then plug it in and switch it on.
3. Wait 20 seconds to allow the router to warm up
4. Switch on the mini routers on the tables
5. Switch on the cameras
6. Run the software in teacher's PC
7. Run the Apps in iPad/Android



Meiji Techno America

5895 Rue Ferrari, San Jose, CA 95138

Phone: 1.800.832.0060 | Fax: 408.226.0900

www.meijitechno.com