
How to upgrade the HD1500MET & HD1500MET-M

camera firmware

Firmware Upgrade steps:


1. Connect the SD card to a PC and create a folder “update” in the SD card root directory. Copy the firmware file “*.bin” to this “update” folder.



Note: The created folder name in the **SD card ROOTDirectory** MUST be “update”.

2. Insert the SD card into the camera SD card slot. Plug in the camera power, attach the USB mouse to the camera. Attach the camera to a monitor via HDMI cable and switch the monitor to HDMI input mode.

<Note: the monitor setting interface might be different because of different brand monitors.>

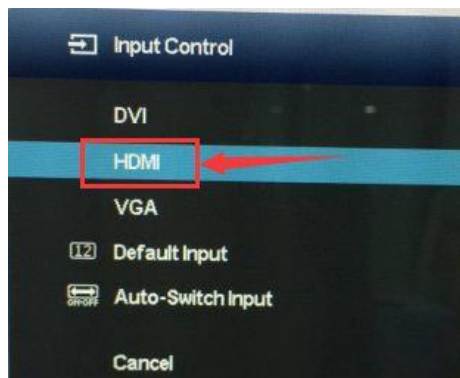

3. Switch on the camera power. A message “Detects new firmware Update” will be popped up in the preview screen, click on “Yes” to start the firmware upgrade.

Note: During the upgrade, please **DO NOT** power off the camera.

4. After finish the upgrade, a message “Update success restart now” shows up. Wait for the camera automatically restarting. Then if it goes to the preview images, it means the firmware is upgraded successfully.

